
AMPTEK Inc.

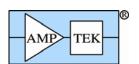


14 DE ANGELO DRIVE, BEDFORD, MA 01730 U.S.A. TEL: (781)-275-2242 FAX: (781)-275-3470 email: sales@amptek.com www.amptek.com


Vacuum Operation

The XR-100 can be operated in air or in vacuum down to 10^{-8} Torr. There are two ways the XR-100 can be operated in vacuum: 1) The entire XR-100 detector and preamplifier box can be placed inside the chamber. In order to avoid overheating and dissipate the 1 Watt of power needed to operate the XR-100, good heat conduction to the chamber walls should be provided by using the four mounting holes. An optional Model 9DVF 9-Pin D vacuum feedthrough connector on a Conflat is available to connect the XR-100 to the PX2 outside the vacuum chamber. 2) The XR-100 can be located outside the vacuum chamber to detect X-Rays inside the chamber through a standard Conflat compression O-ring port. Optional Model EXV9 (9 inch) vacuum detector extender is available for this application. See photograph of XR-100 with extender and Conflat.


Both Option A & B for vacuum chamber pressures up to 10 $^{-8}$ torr OPTION A

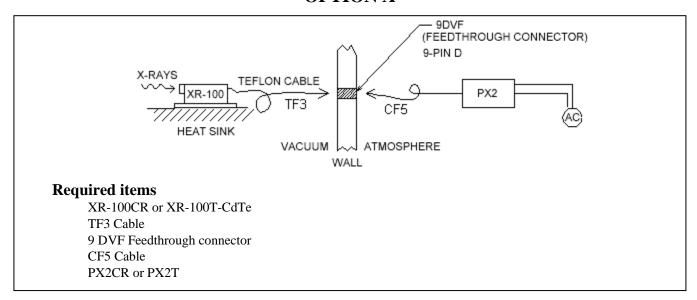

OPTION B

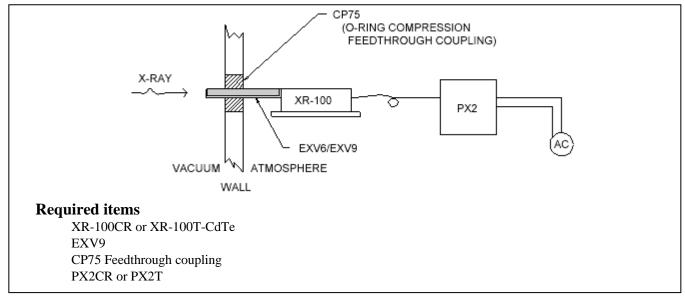

Option B Mechanical Dimensions

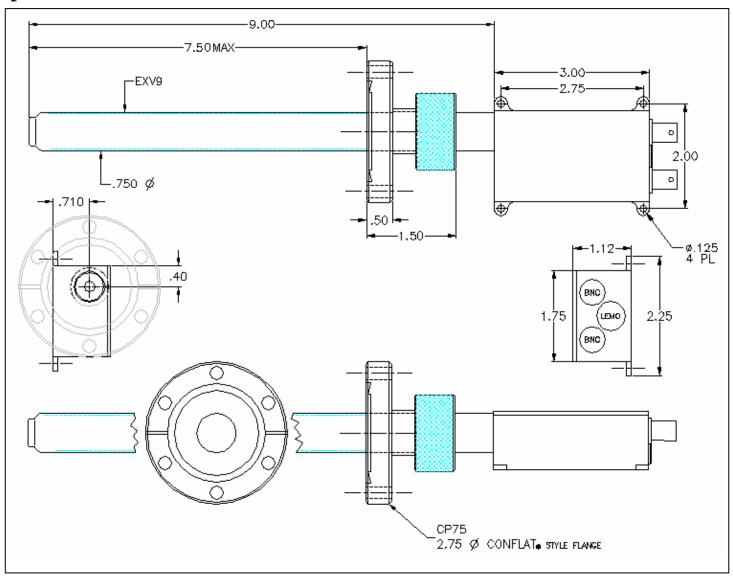





VACUUM APPLICATION ACCESSORIES

Model EXV9


9 Inch Detector Vacuum Extension for the XR-100CR or XR-100T-CdTe

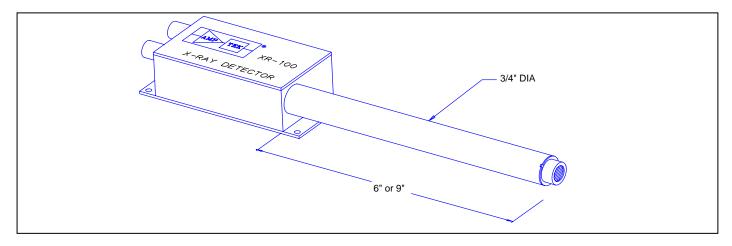


Model CP75
Vacuum Feed Through Coupling on 2 3/4 Inch S.S. Conflat for use with EXV9

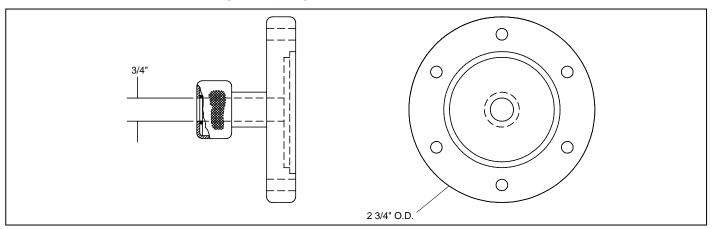

Model 9DVF
9 Pin D-Subminiature Vacuum Feedthrough Connector mounted on a standard 2 3/4 inch S.S. Conflat

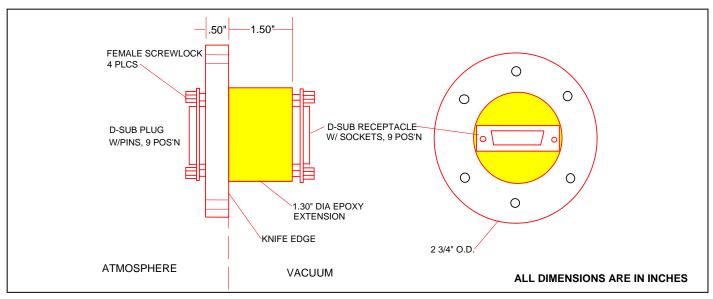

EXVC COLLIMATOR KIT FOR THE EXV9

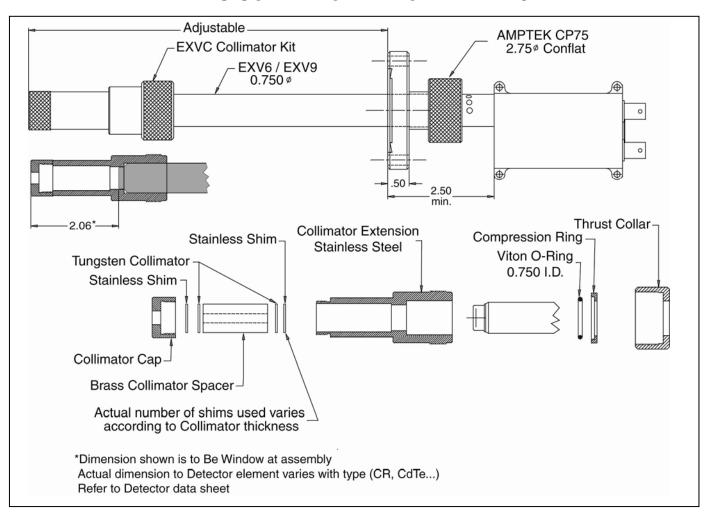

The EXVC - Collimator Kit Includes

Stainless steel collimator housing

Brass Spacer

Tripod and mounting plate (included when ordering for use with standard 1.5 inch extender box) Laser pointer

7 Tungsten (W) Collimator disks:

- 0 1 mm thick with 25 μm hole
- o 1 mm thick with 50 μm hole
- o 2 mm thick with 100 μm hole
- o 2 mm thick with 200 µm hole
- o 2 mm thick with 400 µm hole
- o 2 mm thick with 1000 μm hole
- o 2 mm thick with 2000 μm hole

Optional: EXVC-W-SPACER

This Tungsten (W) Spacer /Collimator is 35 mm thick with a 350 µm hole. It is designed to stop and collimate x-rays greater than 100 keV produced from high energy tubes.

All Tungsten disks are made of alloy HD17 (90% W, 6% Ni, 4% Cu).

